Bedienungsanleitung
Operating Instructions
Mode d'emploi

- » Maxi- und Big-Star Exsikkatoren
- » Maxi- and Big-Star Desiccators
- » Dessiccateurs Maxi- et Big-Star

Sicco Maxi- und Big-Star Exsikkatoren

SICCO Exsikkatoren, Artikelnummer: V 1760-07, V 1835-07, V 1845-07, V 1852-01, V 1853-01, V 1896-07, V 1922-07, V 1923-07, V 1925-07, V 1927-07, V 1844-07, V 1846-07, V 1939-07, V 1936-07, V 1937-07, V 1926-07, V 1928-07 und V 1897-07

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OPERATING INSTRUCTIONS

1. Supplied with	8
2. Security Advice	8
3. Taking into operation	9
4. Functionality	10
5. Service / Maintenance	11
6. Disturbances / Repairs	11
7. Waste disposal	12
8. Specifications	12
9. Spare Parts / Accessories	12

SICCO Big-Star-Desiccators

Versions:

Big-Star-Desiccator
Big-Star-Desiccator Antistatic
Big-Star-Vitrum-Desiccator
Big-Star-Desiccator Black
Big-Star-Desiccator Protect
Big-Star-Desiccator Low

SICCO Maxi 1-Desiccators

Versions:

Maxi 1-Desiccator
Maxi 1-Desiccator Antistatic
Maxi 1-Vitrum-Desiccator
Maxi 1-Desiccator Black
Maxi 1-Desiccator Protect
Maxi 1-Desiccator Horizontal

SICCO Maxi 2-Desiccators

Versions:

Maxi 2-Desiccator
Maxi 2-Desiccator Antistatic
Maxi 2-Vitrum-Desiccator
Maxi 2-Desiccator Black
Maxi 2-Desiccator Protect

SICCO Dolly Desiccator

Sicco Maxi- und Big-Star Desiccators

Please read the information listed below thoroughly. We urgently ask you to observe the instructions about security, use and maintenance.

SICCO Desiccators are made of high-quality plastic material respectively made of transparent Borosilicate glass 3.3. The components of the aluminium frame are joint by connectors made of polyamide.

They are made for safe and secure storage of humidity-sensitive products such as optical and microelectronic parts, reference materials, retained samples or DNA samples in a dry and dust-free atmosphere. This happens by adding SILICA GEL which considerably reduces the air humidity by its absorbing effect.

1. Supplied with

	Shelves made of stainless steel	2 Rails for shelves	Hygrometer	Desiccant tray	Operating Instructions
Big-Star-Desiccators	2	4	1	1	1
Big-Star-Desiccator Low	1	2	1	1	1
Maxi 1-Desiccators	4	8	1	1	1
Maxi 1-Desiccator Horizontal	2	4	1	1	1
Maxi 2-Desiccators	4	8	2	2	1
Dolly-Desiccator	4	8	1	1	1

2. Security Advice


- » Do not take a damaged desiccator into operation.
- » Do not store any solvents or acids in the desiccator.
- » Do not put any explosive or inflammable products into the desiccator.
- » SICCO Desiccators are not suitable for overpressure or vacuum.
- » Provide a safe stand for the desiccator.
- » Place the desiccator in a reachable working height.
- » The ambient temperature range of -20 °C to $+70\text{ °C}$ respectively of -70 °C to $+150\text{ °C}$ for models Vitrum shall not be exceeded.
- » Do not expose the desiccator to direct sunlight.
- » Do not remove more than one shelf at once.

3. Taking into operation

» Remove the protection cover.

» Remove the protection film from the rails and the shelves.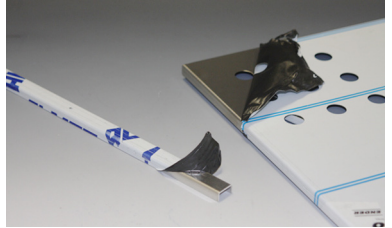


» Opening the door: Hold the desiccator housing with one hand while you open the door by pulling the door grip with the other hand.

» **Please note:** Due to the strong magnetic lock, it is harder to open the door the longer it was closed before. This is normal and can be compensated by pulling stronger than in everyday use.

» Hook the rails (2) into the lateral mounted suspension arrangements (snap-in pin (1) up wards).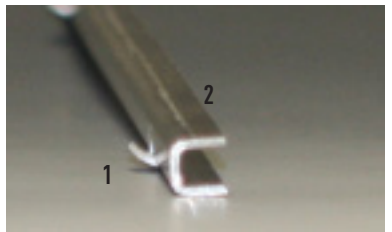

» Mount the rails (2) horizontally in pairs. Pay attention to mount them in the same height. A slight rotation of the rail helps to fix the snap-in pin safely in the suspension arrangement (3).


» For models with rail (2) and suspension arrangement (3) made of plastic: Mount the rails (2) horizontally in pairs on both sides of the suspension arrangement (3). The engagement of the rail (2) is clearly audible. For a better orientation, the levels are marked by numbers on the suspension arrangement (3).

Sicco Maxi- und Big-Star Desiccators


» The shelves can then be slid in on the rails.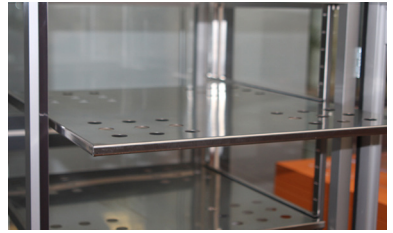

» Fill the tray (available as accessory) with silica gel. The desiccant can be used immediately.


- » Open the battery case on the back of the hygrometer by pushing the cover downwards.
- » Remove the battery safety strip and reinsert the battery. Close the battery case.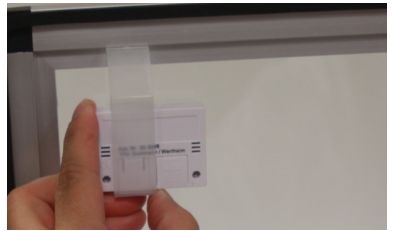

The following function is only available with SICCO Maxi-Desiccators Antistatic and Big-Star-Desiccator Antistatic Cat.-No. V 1922-07, V 1923-07 and V 1925-07.

» Connect the ground cable on the back of the desiccator.

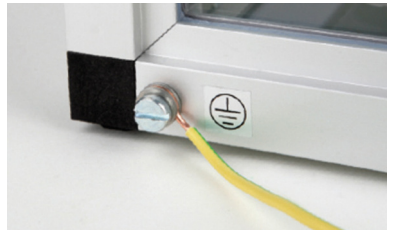4. Functionality

The included silica gel is very hygroscopic and absorbs all air humidity inside the desiccator. The silica gel contains a colour indicator. In unsaturated state, the silica gel is orange. With increasing saturation, the colour is changing from dark brown to black. The silica gel can be regenerated by putting it into an oven for 90 minutes at about 90 to 110 °C. The dumping height influences the time for regeneration: each centimetre in height will take about 90 minutes for regeneration. A microwave oven is not suitable!

Functions of the digital Thermometer/Hygrometer

Measuring range:	<ul style="list-style-type: none">» Temperature – 10 °C to + 60 °C.» Air humidity 20 % to 90 % relative humidity.
------------------	--

Sicco Maxi- und Big-Star Desiccators

Max. error in measurement:	<ul style="list-style-type: none"> » +/- 1°C at -10 °C to + 50 °C. » +/- 5 % relative humidity at 25 % to 75 % relative humidity and 0 °C to + 50 °C. » Attention: The mounted thermometer/hygrometer cannot display any values or data if the relative inside humidity of the desiccator is less than 25 % or higher than 75 %. To display exact values, we recommend using a thermometer/hygrometer of higher quality.
Batterie:	<ul style="list-style-type: none"> » Please use a button cell LR44 or AG13.
Functionality of the Keys:	<p>Indication of maximum / minimum values:</p> <ul style="list-style-type: none"> » By pressing the MAX/MIN-key, the maximum value since the last reset will be displayed. » By pressing the MAX/MIN-key again, the minimum value since the last reset will be displayed. » The current values will be displayed after pressing MAX/MIN a third time. <p>Reset of the MAX/MIN-memory:</p> <ul style="list-style-type: none"> » Press CLEAR and the MAX/MIN-memory will be deleted. The current values will be stored instead. <p>Battery sign in the display</p> <ul style="list-style-type: none"> » If the battery is low, a battery flag is shown on the left bottom of the display. Exchange the battery to ensure best possible operation of the device. <p>SICCO Super-Star-Desiccator (V1875-07) only:</p> <ul style="list-style-type: none"> » The key on the back of the thermometer/hygrometer allows to display the temperature either in °C or °F.

5. Service / Maintenance

- » Use only soft clothes for cleaning. Do not use any detergents which may contain solvents or which scrub.
- » Keep the magnetic door strip clean.

6. Disturbances / Repairs

The following information shall help you to recognise and, if possible, to eliminate possible reasons for disturbances by yourself. If you are not able to handle the disturbance, please give us a call. We will then try to help you this way. If then, a solution to your problem cannot be found, we will ask you to return the cleaned desiccator into the original packing to us. Then we will check, whether a cost-effective repair is possible.

Objection	Reason and help
Humidity inside the desiccator does not fall.	<ul style="list-style-type: none"> » The connecting pieces are not assembled properly. Check again. » The door is not closed tightly. Close the door. » The battery in the hygrometer is empty. Exchange the battery. » The silica gel is saturated. Regenerate the silica gel as described under point 4. » Perhaps you have not added enough silica gel. Add further silica gel. » The built-in hygrometer might be faulty. Double-check the displayed values with an external hygrometer. Exchange the hygrometer.

Sicco Maxi- und Big-Star Desiccators

7. Waste Disposal

The desiccators as well as their packing are made of recyclable materials. Please take further information on the materials used from the following specifications.

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BOHLENDER GmbH